

NucleoSpin[®] Soil

February 2023/Rev. 12



MACHEREY-NAGEL

www.mn-net.com

Contact MN

Germany and international

MACHEREY-NAGEL GmbH & Co. KG Valencienner Str. 11 · 52355 Düren · Germany Tel.: +49 24 21 969-0 Toll-free: 0800 26 16 000 (Germany only) E-mail: info@mn-net.com

Technical Support Bioanalysis

Tel.: +49 24 21 969-333 E-mail: support@mn-net.com

USA

MACHEREY-NAGEL Inc. 924 Marcon Blvd. · Suite 102 · Allentown PA, 18109 · USA Toll-free: 888 321 6224 (MACH) E-mail: sales-us@mn-net.com

France

 MACHEREY-NAGEL SAS

 1, rue Gutenberg – BP135 · 67720 Hoerdt Cedex · France

 Tel.:
 +33 388 68 22 68

 E-mail:
 sales-fr@mn-net.com

 MACHEREY-NAGEL SAS (Société par Actions Simplifiée) au capital de 186600 €

 Siret 379 859 531 0020 · RCS Strasbourg B379859531 · N° intracommunautaire FR04 379 859 531

Switzerland

MACHEREY-NAGEL AG Hirsackerstr. 7 · 4702 Oensingen · Switzerland Tel.: +41 62 388 55 00 E-mail: sales-ch@mn-net.com

www.mn-net.com

Genomic DNA from soil

Protocol at a glance (Rev. 12)

NucleoSpin® Soil

| 1 | Prepare sample | MN Bead Tube Type A | | | | |
|----|---------------------------|---------------------|--|--|--|--|
| | | | | 250 – 500 mg sample material | | |
| | | M | | 700 µL SL1 or SL2 | | |
| 2 | Adjust lysis conditions | | 150 µL Enhancer SX | | | |
| 3 | Sample lysis | | | Horizontally vortex 5 min at RT or use other homogenizers according to manufacturers protocol | | |
| 4 | Precipitate | | | 11,000 × <i>g</i> , 2 min | | |
| | contaminants | | | 150 µL SL3 | | |
| | | | Ò | Vortex 5 s | | |
| | | M | | 0-4 °C, 5 min | | |
| | | | | 11,000 × <i>g</i> , 1 min | | |
| 5 | Filter lysate | | Ċ | Load supernatant on NucleoSpin® Inhibitor Removal Column (red ring) | | |
| | | | | 11,000 × <i>g</i> , 1 min | | |
| 6 | Adjust binding conditions | | | 250 µL SB | | |
| | | U | | Vortex 5 s | | |
| 7 | Bind DNA | Ĵ | Load 550 µL sample NucleoSpin® Soil Column (c | | | |
| | | | Ć | 11,000 × <i>g</i> , 1 min | | |
| | | | 0 | Load remaining sample | | |
| | | | | 11,000 × <i>g</i> , 1 min | | |
| 8 | Wash silica membrane | | | 1 st 500 μL SB 11,000 × <i>g</i> , 30 s | | |
| | membrane | | | 2nd 550 μL SW1 11,000 × <i>g</i> , 30 s | | |
| | | | \bigcirc | 3 rd 650 µL SW2 Vortex 2 s 11,000 x <i>g</i> , 30 s | | |
| | | | | 4 th 650 μL SW2 Vortex 2 s 11,000 x <i>g</i> , 30 s | | |
| 9 | Dry silica | Ĩ | | | | |
| | membrane | | Ò | 11,000 x <i>g</i> , 2 min | | |
| 10 | Elute DNA | | | 30–100 µL SE | | |
| | | | Ć | RT, 1 min | | |
| | | ¥ | Ű | 11,000 x <i>g</i> , 30 s | | |
| | | | | , | | |



Table of contents

| 1 | Con | nponents | 4 |
|---|------|---|----|
| | 1.1 | Kit contents | 4 |
| | 1.2 | Reagents, consumables, and equipment to be supplied by user | 5 |
| | 1.3 | About this user manual | 5 |
| 2 | Pro | duct description | 6 |
| | 2.1 | The basic principle | 6 |
| | 2.2 | Kit specifications | 6 |
| | 2.3 | Relevance of humic substances as PCR inhibitors | 7 |
| | 2.4 | Amount of starting material | 8 |
| | 2.5 | Choice of lysis buffer | 8 |
| | 2.6 | Mechanical sample lysis | 10 |
| | 2.7 | Repeated extraction | 10 |
| | 2.8 | Elution procedures | 10 |
| | 2.9 | How to interpret DNA yield and purity from UV-VIS | 11 |
| | 2.10 | Isolation of DNA from water samples | 13 |
| | 2.11 | Isolation of DNA from samples stabilized in Zymo DNA/RNA Shield | 15 |
| 3 | Stor | age conditions and preparation of working solutions | 16 |
| 4 | Safe | ety instructions | 17 |
| | 4.1 | Disposal | 17 |
| 5 | Prot | ocol – purification of DNA from soil and sediment | 18 |
| | 5.1 | Protocol – purification of DNA from water | 22 |
| 6 | Арр | endix | 24 |
| | 6.1 | Troubleshooting | 24 |
| | 6.2 | Ordering information | 26 |
| | 6.3 | Product use restriction/warranty | 28 |

1 Components

1.1 Kit contents

| | NucleoSpin [®] Soil | | |
|--|------------------------------|-----------------------|-------------------------|
| REF | 10 preps 740780.10 | 50 preps 740780.50 | 250 preps 740780.250 |
| Lysis Buffer SL1 | 30 mL | 60 mL | 250 mL |
| Lysis Buffer SL2 | 30 mL | 60 mL | 250 mL |
| Lysis Buffer SL3 | 10 mL | 10 mL | 50 mL |
| Enhancer SX | 3 mL | 10 mL | 50 mL |
| Binding Buffer SB | 10 mL | 60 mL | 250 mL |
| Wash Buffer SW1 | 6 mL | 30 mL | 150 mL |
| Wash Buffer SW2 (Concentrate)* | 6 mL | 25 mL | 100 mL |
| Elution Buffer SE** | 13 mL | 13 mL | 60 mL |
| MN Bead Tubes Type A | 10 | 50 | 250 |
| NucleoSpin [®] Inhibitor Removal Columns (red rings) | 10 | 50 | 250 |
| NucleoSpin [®] Soil Columns (green rings) | 10 | 50 | 250 |
| Collection Tubes (2 mL) | 10 | 50 | 250 |
| Collection Tubes (2 mL, lid) | 10 | 50 | 250 |
| User manual | 1 | 1 | 1 |

^{*} For preparation of working solutions and storage conditions see section 3.

^{**} Composition of Elution Buffer SE: 5 mM Tris/HCl, pH 8.5

1.2 Reagents, consumables, and equipment to be supplied by user

Reagents

- Isopropanol (when working with samples which are stabilized in Zymo DNA/RNA Shield)
- 96-100 % ethanol

Consumables

- 1.5 mL microcentrifuge tubes
- Disposable pipette tips

Equipment

- Manual pipettors
- Centrifuge for microcentrifuge tubes
- Equipment for sample disruption and homogenization (see section 2.6)
- Personal protection equipment (e.g., lab coat, gloves, goggles)

1.3 About this user manual

It is strongly recommended that first time users of the **NucleoSpin[®] Soil** kit read the detailed protocol sections of this user manual. Experienced users, however, may refer to the Protocol at a glance instead. The Protocol at a glance is designed to be used only as a supplemental tool for quick referencing while performing the purification procedure.

All technical literature is available on the Internet at *www.mn-net.com*.

Please contact Technical Service regarding information about changes of the current user manual compared to previous revisions.

2 Product description

2.1 The basic principle

The sample material is resuspended in Lysis Buffer SL1 or SL2, supplemented with the Enhancer SX, and mechanically disrupted using ceramic beads.

Proteins and PCR inhibitors are precipitated with Lysis Buffer SL3 and subsequently pelleted by centrifugation together with the ceramic beads and undissolved sample material. The supernatant is taken off and cleared by passing it through a NucleoSpin[®] Inhibitor Removal Column.

DNA binding conditions are then adjusted by addition of Binding Buffer SB to the flow through and the lysate is loaded onto a NucleoSpin[®] Soil Column.

Residual humic substances, especially humic acids, and other PCR inhibitors are removed by efficient washing with Binding Buffer SB and Wash Buffers SW1/SW2. After a drying step, ready to use DNA can be eluted with Elution Buffer SE (5 mM Tris/HCl, pH 8.5).

2.2 Kit specifications

- The NucleoSpin[®] Soil kit is designed for the isolation of high molecular weight genomic DNA from microorganisms like Gram positive and Gram negative bacteria, archaea, fungi, and algae in soil, sludge, and sediment samples.
- Suitable for soils from forest, bog, farmland, grassland, etc.
- Suitable for samples stabilized in Zymo DNA/RNA Shield (isopropanol required)
- Suitable for stool samples.
- The kit offers two special lysis buffers, **Buffer SL1** and **Buffer SL2**, which can be combined with the chemical additive **Enhancer SX** to guarantee highest possible yields with excellent purity for all types of sample material.
- Efficient mechanical lysis of the sample material is achieved by bead beating using the ceramic **MN Bead Tubes Type A**.
- The optimized buffer chemistry and the NucleoSpin[®] Inhibitor Removal Column completely remove humic substances and other PCR inhibitors typically present in soil and sediment samples.
- The eluted DNA is ready to use for all standard downstream applications. In most cases the concentrated DNA can be used as PCR template without further dilution for highest sensitivity.

| Table 1: Kit specifications at a glance | | | |
|---|------------------------------|--|--|
| Parameter | NucleoSpin [®] Soil | | |
| Format | Mini spin columns | | |
| Sample material | < 500 mg soil or sediment | | |
| Typical yield | 2–10 µg | | |
| Elution volume | 30–100 μL | | |
| Preparation time | 90 min/10 preps | | |
| Binding capacity | 50 µg | | |
| Use | For research use only | | |

2.3 Relevance of humic substances as PCR inhibitors

Humic substances are produced by bacteria, fungi, and protozoa in soil, sediments and waters during the degradation of plant or other organic matter. They consist of very high molecular weight compounds with undefined structures. Building blocks are mainly heterocyclic aromatic compounds that are linked by ether or ethoxy groups and which carry hydroxyl-, methoxy-, carbonyl-, or carboxyl groups.

According to their solubility in water they are divided into humin, humic acids, and fulvic acids. The completely insoluble and black humin has an average molecular weight of around 300,000 g/mol. The dark brown to grey colored humic acids are slightly smaller. They carry a lot of hydroxyl and carboxyl groups and are therefore mainly soluble at neutral or alkaline pH. The only slightly yellow to light-brown colored fulvic acids with an average molecular weight of 2,000 g/mol are soluble under alkaline as well as under acidic conditions.

Due to the high molecular weight and the mainly polyanionic nature of humic substances most purification methods do not distinguish between these molecules and DNA. For the same reason they act as extremely potent PCR inhibitors. Even smallest amounts of humic substances can inhibit for example DNA polymerases or restriction enzymes and result in a complete failure of enzymatic downstream applications.

Frequently, the problem is circumvented by dilution of the isolated DNA prior to PCR analysis. However, this results in a significantly reduced sensitivity because low abundance DNA may be lost completely.

Thus, highest DNA yields with as little PCR inhibitor contaminations as possible are of utmost importance for any DNA analysis of soil samples.

2.4 Amount of starting material

NucleoSpin[®] Soil is suitable for processing 250 – 500 mg of sample material. However, do not fill the MN Bead Tube Type A higher than the 1 mL mark (including the ceramic beads) to ensure sufficient head space for an efficient mechanical disruption.

Usually a reduction of starting material also helps to improve the lysis efficiency and to increase the purity of the DNA.

Very dry material can soak up large volumes of lysis buffer. In this case, either reduce the amount of sample material or add additional lysis buffer up to the 1.5 mL mark of the MN Bead Tube Type A.

If possible remove foreign material like leaves, stones, or twigs (e.g., by sieving) as well as excess of water (e.g., by discarding the supernatant after spinning down sediment samples).

2.5 Choice of lysis buffer

Due to the highly varying composition of different soils (organic matter, inorganic matter, humic substances, metal ions, polysaccharides, pH, etc.), it is impossible to obtain best results in DNA yield and purity for all sample types with only one single lysis buffer system.

There are several parameters that can be adjusted in a way that lysis works perfect for one sample but fails with another. Therefore, the NucleoSpin[®] Soil kit is equipped with two lysis buffers SL1 and SL2 and an Enhancer SX.

Those three components allow a perfect fine tuning for every type of soil sample for maximum yield and purity. Unfortunately, for the reasons given above there is no way to predict the best choice of lysis buffer for a specific sample. This can only be determined experimentally. Therefore, **both lysis buffers should be tested in parallel** for each new sample material. Usually Buffer SL1 in combination with Enhancer SX is more suitable for soil consisting predominantly of minerals while Buffer SL2 is more suitable for soil with a high amount of organic carbon.

After mixing the sample with lysis buffer in the MN Bead Tube Type A, the Enhancer SX is added routinely to the sample prior to the mechanical homogenization. This buffer ensures the highest possible DNA yield with most sample materials. However, in case of a very high humic acid content in the sample material, the Enhancer SX might also reduce the purity of the DNA by facilitating the release of humic acids into the lysate. Therefore, the volume of added Enhancer SX can be lowered from 150 μ L to for example 10 μ L or the buffer can be entirely omitted. This usually increases the purity (A₂₆₀/A₂₃₀) of the sample significantly (Table 2), might, however, lower the DNA yield (Figure 1).

Ideally, for a new sample material both lysis buffers **Buffer SL1 and SL2** should be tested **with and without adding Enhancer SX.** These initial four preparations will help you to find the ideal lysis condition for your special soil composition.

| | - | | | |
|----------------|---|---|---|---|
| 1 | 2 | 3 | 4 | 5 |
| - | _ | _ | _ | - |
| | | | | |
| <u>heat-of</u> | | | | |
| | | | | |

Figure 1:Total DNA purified from wheat field soil with four different lysis buffer combinations

20 of 100 μ L eluate were analyzed on a 1 % TAE agarose gel:

Lane 1: Marker λ / HindIII

Lane 2: Lysis Buffer SL1

Lane 3: Lysis Buffer SL1 +Enhancer SX

Lane 4: Lysis Buffer SL2

Lane 5: Lysis Buffer SL2 +Enhancer SX

| Table 1: Yields and purity ratios of DNA purified from wheat field soil | | | | | |
|---|--------|--------|--------|--------|--|
| Buffer | SI | L1 | SL2 | | |
| Enhancer SX | - | + | - | + | |
| Yield | 2.3 µg | 2.3 µg | 1.4 µg | 3.1 µg | |
| A ₂₆₀ /A ₂₈₀ | 1.69 | 1.60 | 1.76 | 1.72 | |
| A ₂₆₀ /A ₂₃₀ | 1.85 | 0.96 | 1.78 | 0.99 | |

2.6 Mechanical sample lysis

A thorough mechanical lysis step is essential to break up the soil crumbs, to free the cells within the soil, and to break up cells and spores. Ceramic beads have proven to be most effective in combination with a bead mill, a FastPrep[®]-24 instrument (MP Biomedicals, set instrument to 5 m/s for 30 s), or an MN Bead Tube Holder, see ordering information. In most cases, however, this kind of equipment is not necessary. The same result can be achieved by taping the lysis tubes **horizontally** to a standard vortexer.

The lysis time should be as short as necessary to avoid shearing of DNA and to minimize the release of humic acids. Depending on the sample, however, it might be advantageous to increase the lysis time to 10, 20, or 30 min.

Homogenization and cell disruption should be performed at room temperature $(18-25 \ ^{\circ}C)$ to avoid SDS precipitation in the lysis buffers. Overheating the sample, for example by prolonged bead beating in a bead mill or the FastPrep[®]-24 instrument, should be avoided to minimize liberation of humic acids.

2.7 Repeated extraction

For sample materials containing a high amount of microorganisms a single extraction step might not be sufficient to disrupt every cell and to release all DNA. Extracting the sample twice may help to increase DNA yield significantly.

Therefore, follow the protocol until the first centrifugation in step 4. But instead of adding SL3 directly to the MN Bead Tube Type A, transfer the supernatant to a new collection tube (not provided) and complete step 4 with this supernatant. Then repeat steps 1-4 with the same soil sample in the MN Bead Tube Type A. Filter both final supernatants of step 4 through a NucleoSpin[®] Inhibitor Removal Column as described in step 5. Add Binding Buffer SB to both filtrates according to step 6 and finally load both samples on one NucleoSpin[®] Soil Column according to step 7 in multiple loading steps.

Note that the supplied buffer volumes are calculated for only one extraction. The excess of Enhancer SX and Binding Buffer SB might not be sufficient to allow two extraction steps for all 10, 50, or 250 preps of the kit.

For more information, please visit our website: https://www.mn-net.com/de/microbiome-information

2.8 Elution procedures


It is possible to adapt the elution method, temperature, and volume of elution buffer used for the subsequent application of interest. In addition to the standard method where an increase of DNA concentration can be achieved by reducing the elution volume from 100 to 30 μ L, there are two options to increase the DNA yield:


- Heat the elution buffer to 80 °C.
- Perform two subsequent elution steps with fresh elution buffer.

2.9 How to interpret DNA yield and purity from UV-VIS

The most common method to determine the DNA yield is UV-VIS spectroscopy. The DNA concentration in the final eluate can be calculated from its absorption maximum at 260 nm (A₂₆₀) based on the fact that an absorption of A₂₆₀ = 1 corresponds to 50 µg/mL double stranded DNA. However, this calculation assumes the absence of any other compound that absorbs UV light at 260 nm. Any contamination with, for example, RNA, protein, or especially humic substances significantly contributes to the total absorption at 260 nm and therefore leads to an overestimation of the real DNA concentration.

Figure 2 shows a typical UV absorbance spectrum of pure DNA (solid line) exhibiting a peak at 260 nm, a decrease of absorption with a minimum at 230 nm, and only a moderate increase in absorption below 230 nm. In comparison, the spectrum of a sample that is contaminated with humic acids demonstrates only a small shoulder at 260 nm, it lacks the minimum at 230 nm, and the absorption sores up below 230 nm. In this case only a small part of the absorbance at 260 nm is caused by DNA, most of it is just the tailing absorption of the humic acid contamination. However, the calculated DNA yield seems to be higher in the contaminated sample. Thus, DNA yield determined by UV-VIS, might be distorted by co-purifying contaminants and we recommend to check the DNA yield also by agarose gel electrophoresis.

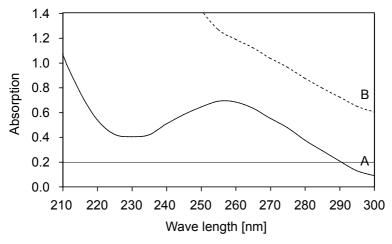

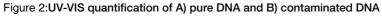

A) 7.7 μg in 100 $\mu L,$ 1.84 $A_{260}/A_{280},$ 1.71 A_{260}/A_{230}

B) 9.3 μg in 100 $\mu L,$ 1.35 $A_{260}/A_{280},$ 0.27 A_{260}/A_{230}

Purity ratio A₂₆₀/A₂₃₀

To facilitate the decision whether the yield as determined from A_{260} readings can be trusted or not, the ratio of the absorption at 260 nm and 230 nm can be used. The ratio A_{260}/A_{230} should be higher than 2.0 for pure DNA and is acceptable down to ratios of about 1.5. Smaller values around or even below 1.0, as shown in Figure 2, indicate significant amounts of impurities and the real DNA concentration is far below its calculated value.

Additionally, not only humic acids, but also proteins, saccharides, and other contaminants can be detected by a low A_{260}/A_{230} ratio.

Purity ratio A₂₆₀/A₂₈₀

Another indicator of DNA purity is the ratio A_{260}/A_{280} , which should be between 1.8 and 1.9. Values below 1.8 indicate protein contamination, whereas higher values indicate RNA contamination. However, this ratio should be treated with caution, since contamination with protein and RNA at the same time can compensate each other and result in a perfect A_{260}/A_{280} .

Agarose gel electrophoresis

As a consequence, the DNA should always be run on an agarose gel to verify the UV-VIS quantification especially if A_{260}/A_{230} and A_{260}/A_{280} are beyond the acceptable range. Figure 3 demonstrates that the contaminated sample B) of Figure 2 actually contains much less DNA than the pure sample A) in contrast to the UV-VIS results, which can easily be misinterpreted.

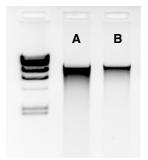


Figure 3:Gel analysis of A) pure and B) contaminated genomic DNA from soil

10 μ L of each sample were run on a 1 % TAE agarose gel (1 h, 100 V). The larger gel band of pure DNA A) proves a higher yield and concentration compared to the contaminated DNA sample which is in contrast to the UV-VIS quantification (A: 7.7 μ g/100 μ L, B: 9.3 μ g/100 μ L).

2.10 Isolation of DNA from water samples

Additionally required equipment

- MN Bead Tubes 5 mL Type A for 45-50 mm filters (see ordering information)
- MN Bead Tube Holder 5 mL (see ordering information)
- 15 mL centrifuge tubes (for use with 45 50 mm filters)

Volume of water sample

The maximum volume of water sample that can be processed varies depending on the sample (e.g. source and quality) and filter membrane (e.g., type, diameter, pore size).

The turbidity of water samples can vary from clear to highly turbid, due to different concentrations of particulates. In general, high volumes of clear or potable drinking water can be processed, as there is a lower chance of filter clogging. A volume of 100 mL up to several liters might be processed.

Turbid water samples containing high levels of sediments or suspended particles, such as clay, silt, or other inorganic or organic matter, may lead to filter clogging. The use of a 0.45 μ m filter is recommended for these samples types. Furthermore, it is advantageous to let samples sediment over time. In case a sedimentation is not desired or does not occur, a pre-filter on top of the membrane filter is recommended. MACHEREY-NAGEL offers a broad range of filter paper suitable for pre-filtration of large debris.

| Suitable pre- filter | Properties | Particle retention | Diameter | REF |
|-------------------------|-------------------|--------------------|----------|--------|
| MN 619 G | slow filtration | approx. 2–7 µm | 55 mm | 440005 |
| MN 616 G | medium filtration | approx. 4–12 µm | 55 mm | 483005 |
| MN 617 G | fast filtration | approx. 7–12 µm | 55 mm | 494005 |

Storage of water sample

The most convenient way is the storage of filtered water samples. Therefore, filter the water sample through a membrane filter and store the filter directly within the appropriate amount of lysis buffer in the respective bead tube. Keep in mind that sodium dodecylsulfate in Buffer SL1 and SL2 will precipitate at lower temperatures. Incubate tubes at 37 °C until the SDS precipitates have redissolved before starting the preparation.

Filter Membranes

NucleoSpin[®] Soil is suitable for the nucleic acid extraction from both 25 mm and 47 mm filter membranes of different composition or origin. Disposable filter funnel units as well as reusable filter funnels are compatible and can be used with different filter membrane types. MACHEREY-NAGEL offers a broad range of different membrane filter types and sizes.

| Туре | Diameter | Pore size | Product | REF |
|------------------------|-----------------|------------------|-------------------------|-------------|
| Cellulose acetate (CA) | 25 mm | 0 . 2 µm | PORAFIL [®] CA | 68000020025 |
| | | 0 . 45 µm | PORAFIL [®] CA | 68000045025 |
| | 47 mm | 0 . 2 µm | PORAFIL [®] CA | 68000020047 |
| | | 0 . 45 µm | PORAFIL [®] CA | 68000045047 |
| Cellulose mixed esters | 25 mm | 0 . 45 µm | PORAFIL [®] CM | 65100045025 |
| (CM) | 47 mm (sterile) | 0 . 45 µm | PORAFIL [®] CM | 65300045047 |
| Cellulose nitrate (NC) | 25 mm | 0 . 2 µm | PORAFIL [®] NC | 6570020025 |
| | | 0 . 45 µm | PORAFIL [®] NC | 6570045025 |
| | 47 mm | 0 . 2 µm | PORAFIL [®] NC | 6570020047 |
| | | 0 . 45 µm | PORAFIL [®] NC | 6570045047 |
| Polycarbonate (PC) | 25 mm | 0 . 40 µm | PORAFIL [®] PC | 676040025 |
| Polyethersulfone (PES) | - | - | | - |

The choice of a suitable bead tube type highly depends on the filter membrane size.

- 25 mm filter membranes: MN Bead Tube Type A (0.6-0.8 mm ceramic beads, supplied)
- 47 mm filter membranes: MN Bead Tube 5 mL Type A (0.6–0.8 mm ceramic beads, see ordering information)

2.11 Isolation of DNA from samples stabilized in Zymo DNA/ RNA Shield

The NucleoSpin[®] Soil kit can be combined with Zymo DNA/RNA Shield solution. Add Zymo DNA/RNA Shield to the soil sample in Bead Tube Type A until the sample is barely but completely submerged. Use as little Zymo DNA/RNA Shield solution as possible to keep the total volume low. 500 mg of soil sample usually require the addition of 500 μ L Zymo DNA/RNA Shield.

Follow the protocol to step 6. Adjust binding conditions by addition of 250 μ L Buffer SB according to step 6 and determine the total sample volume. Add 0.2 sample volumes of isopropanol to the sample and vortex for 5 s. For example: if the total sample volume after addition of Buffer SB is 1000 μ L, add 200 μ L isopropanol. Continue with step 7.

3 Storage conditions and preparation of working solutions

Attention:

Buffers SB and SW1 contain guanidinium thiocyanate and guanidine hydrochloride, respectively. Wear gloves and goggles!

Storage conditions:

 All kit components should be stored at 15-25 °C and are stable until: see package label. Storage at lower temperatures may cause precipitation of salts. If precipitation occurs, incubate the bottle for several minutes at about 30-40 °C and mix well until the precipitate is dissolved.

Before starting the first NucleoSpin[®] Soil procedure prepare the following:

- Wash Buffer SW2: Add the indicated volume of ethanol (96–100%) to Buffer SW2 Concentrate. Mark the label of the bottle to indicate that ethanol was added. Buffer SW2 is stable at 15–25 °C for at least one year.
- Sodium dodecyl sulfate (SDS) in Buffer SL1/SL2 may precipitate if stored at temperatures below 20 °C. If a precipitate is observed in Buffer SL1/SL2, incubate the bottle at 30–40 °C for several minutes and mix well

| NucleoSpin [®] Soil | | | | | |
|----------------------------------|---------------------------|--------------------------------|---------------------------------|--|--|
| REF | 10 preps 740780.10 | 50 preps 740780.50 | 250 preps 740780.250 | | |
| Wash Buffer SW2 (Concentrate) | 6 mL Add 24 mL ethanol | 25 mL Add 100 mL ethanol | 100 mL Add 400 mL ethanol | | |

4 Safety instructions

When working with the **NucleoSpin[®] Soil** kit wear suitable protective clothing (e.g., lab coat, disposable gloves, and protective goggles). For more information consult the appropriate Material Safety Data Sheets (MSDS available online at **www.mn-net.com/msds**).



Caution: Guanidin thiocyanate in Buffer SB and buffer SW1 can form highly reactive compounds when combined with bleach! Thus, do not add bleach or acidic solutions directly to the sample preparation waste.

The waste generated with the **NucleoSpin[®] Soil** kit has not been tested for residual infectious material. A contamination of the liquid waste with residual infectious material is highly unlikely due to strong denaturing lysis but it cannot be excluded completely. Therefore, liquid waste must be considered infectious and should be handled and discarded according local safety regulations.

4.1 Disposal

Dispose hazardous, infectious or biologically contaminated materials in a safe and acceptable manner and in accordance with all local and regulatory requirements.

5 Protocol – purification of DNA from soil and sediment

Before starting the preparation:

• Check Lysis Buffer SL1 or SL2 for precipitated SDS. Dissolve any precipitate by incubating the buffer at 30–40 °C for 10 min and shaking the bottle every 2 min.

1 Prepare sample

See section 2.4 and 2.5 for more information on the amount of starting material and the choice of lysis buffer. See section 2.7 for the repeated extraction of a sample to improve DNA yield.

sample +700 µL SL1 or SL2

250 – 500 mg

Important: Do not fill the tube higher than the 1 mL mark.

Transfer 250–500 mg fresh sample material to a MN Bead Tube Type A containing the ceramic beads.

Add 700 µL Buffer SL1 or Buffer SL2.

<u>Note for very dry material:</u> If the sample material soaks up too much lysis buffer, fill the MN Bead Tube Type A up to the 1.5 mL mark with fresh lysis buffer.

<u>Note for very wet material:</u> Remove excess liquid before addition of lysis buffer, if necessary after spinning down the sample.

2 Adjust lysis conditions

Add 150 µL Enhancer SX and close the cap.

<u>Note:</u> Enhancer SX ensures the highest possible DNA yield. It can, however, also promote the release of humic acids. See section 2.5 on how to lower the volume or omit the buffer entirely in order to increase DNA purity.

3 Sample lysis

See section 2.6 for more information on homogenization methods (e.g., FastPrep[®]-24 instrument, Vortex adapter).

Attach the MN Bead Tubes **horizontally** to a vortexer, for example, by taping or using a special adapter.

Vortex the samples at **full speed** and **room temperature** (18-25 °C) for **5 min**.

+150 μL SX

Vortex RT, 5 min

4 Precipitate contaminants

Centrifuge for **2 min** at **11,000 x** *g* to eliminate the foam caused by the detergent.

<u>Note:</u> The clear supernatant can be transferred to a new collection tube (not provided) prior to the following precipitation. This might result in more consistent yields from prep to prep and is highly recommended for carbonate containing samples. See also section 2.7 for repeated extraction of a sample to improve DNA yield.

Add 150 µL Buffer SL3 and vortex for 5 s.

Incubate for 5 min at 0-4 °C.

Centrifuge for **1 min** at **11,000 x** g.

5 Filter lysate

Place a **NucleoSpin[®] Inhibitor Removal Column** (red ring) in a Collection Tube (2 mL, lid).

Load up to $700\;\mu L$ clear supernatant of step 4 onto the filter.

Centrifuge for **1 min** at **11,000 x** *g*.

<u>Note:</u> With very wet samples (e.g., sediments) the volume of clear supernatant of step 4 can exceed 700 μ L significantly. In this case transfer the NucleoSpin[®] Inhibitor Removal Column to a new collection tube (not provided) and load the remaining supernatant. Centrifuge for 1 min at 11,000 x g. Combine the flow throughs.

Discard the NucleoSpin® Inhibitor Removal Column.

If a pellet is visible in the flow through, transfer the clear supernatant to a new collection tube (not provided).

6 Adjust binding conditions

Add 250 µL Buffer SB and close the lid.

Vortex for 5 s.

<u>Note:</u> If samples were stabilized in Zymo DNA/RNA Shield, quantify total sample volume after addition of Buffer SB and add 0.2 volumes of isopropanol. 11,000 x g, 2 min +150 μL SL3 Vortex 5 s 0-4 °C.

5 min



Load supernatant

> +250 μL SB Vortex 5 s

7 Bind DNA

Place a NucleoSpin® Soil Column (green ring) in a Collection Tube (2 mL).

Load 550 µL sample onto the column.


Centrifuge for 1 min at 11,000 x g.

Discard flow through and place the column back into the collection tube.

Load the remaining sample onto the column.

Centrifuge for 1 min at 11,000 x g.

Discard flow through and place the column back into the collection tube.

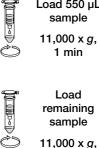


Load remaining sample

Load 550 µL

1 min

11,000 x g, 1 min

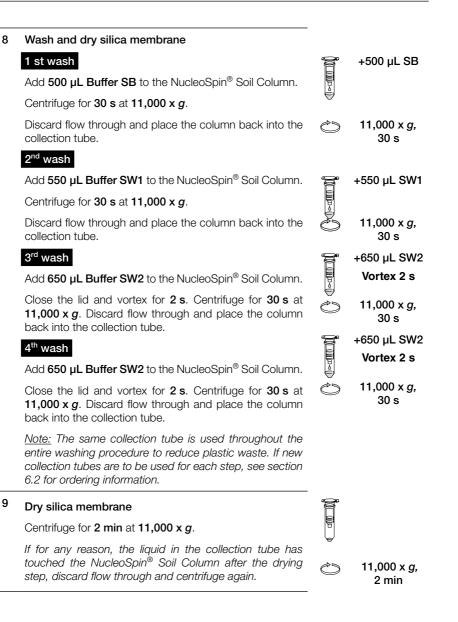

10 Elute DNA

Place the NucleoSpin[®] Soil Column into a new microcentrifuge tube (not provided).

Add $30 \ \mu L$ (for high concentration), $50 \ \mu L$ (for medium concentration and yield), or $100 \ \mu L$ (for high yield) **Buffer SE** to the column.

Do not close the lid and incubate for 1 min at room temperature (18–25 °C). Close the lid and centrifuge for **30 s** at **11,000 x** g.

<u>Note:</u> Quantify DNA not only by UV-VIS but also run an agarose gel to verify yield and DNA quality (see section 2.9 for more information).





5.1 Protocol – purification of DNA from water

Protocol for isolation of DNA from 25 mm Filters

Filter water sample through an appropriate 25 mm filter membrane (not provided, see ordering information) using a suitable filtration device.

Remove the filter membrane from the filtration device by using sterile forceps, roll the membrane into a cylinder (top side facing inwards) and insert into a MN Bead Tube Type A. Add 700 μ L Buffer SL1 or SL2 and optionally up to 150 μ L Buffer SX. Addition of Buffer SX is not recommended for water samples with a high concentration of humic acids.

Continue with step 3 of the protocol. Remove and discard filter using a sterile pipette tip.

Protocol for isolation of DNA from 47 mm Filters

Filter water sample through an appropriate 47 mm filter membrane (not provided, see ordering information) using a suitable filtration device.

Remove the filter membrane from the filtration device by using sterile forceps, roll the membrane into a cylinder (top side facing inwards) and insert it into a MN Bead Tube 5 mL Type A (see ordering information). Add 1400 μ L Buffer SL1 or SL2 and optionally up to 300 μ L Buffer SX. Addition of Buffer SX is not recommended for water samples with a high concentration of humic acids.

Vortex the MN Bead Tube 5 mL Type A on the MN Bead Tube Holder 5 mL (see ordering information) horizontally in conjunction with a Vortex-Genie[®] for 15 min at maximum speed.

Centrifuge the Bead Tubes at > 1000 x g for 2 minutes. Remove and discard filter using a sterile pipette tip.

Add 300 μ L Buffer SL3 and vortex for 5 s. Incubate at 0–4 °C for 5 min.

Centrifuge the Bead Tubes at > 1000 x g for 5 minutes at 4 °C.

Load up to 650 μ L supernatant into a NucleoSpin[®] Inhibitor Removal Column (red ring) and centrifuge at 11,000 x g for 1 min. Collect flow-through in a clean 15 mL centrifuge tube (not supplied). Repeat this step until the remaining supernatant is filtered using fresh NucleoSpin[®] Inhibitor Removal Columns (not supplied, see ordering information) for each repetition.

Add 500 μL Buffer SB to the pooled flow-through and vortex for 5 s. See ordering information for additional Buffer SB if needed.

Load up to 550 μ L sample into a NucleoSpin[®] Soil Column (green ring) and centrifuge at 11,000 x g for 1 min. Discard flow-through. Repeat this step to load the remaining sample.

Continue with step 8 of the protocol.

6 Appendix

6.1 Troubleshooting

| Problem | Possible cause and suggestions |
|-------------------------|---|
| | Suboptimal lysis conditions |
| | • Too much sample material was filled into the MN Bead Tube Type A. Too little head space does not allow the necessary motion of the beads to disrupt the sample. Use less sample material (see section 2.4 for more information). |
| | • Compare the yields obtained with Lysis Buffer SL1 and SL2 in parallel purifications each with and without addition of Enhancer SX to find the optimal lysis buffer conditions (see section 2.5 for more information). |
| | Insufficient disruption and/or homogenization of starting material |
| | • Shaking of the MN Bead Tube Type A was too weak or not long enough. Increase shaking time and velocity or use another shaking device (see section 2.6 for more information). Make sure that the MN Bead Tube Type A is fixed horizontally on the vortexer. |
| | Reagents not applied or restored properly |
| Poor or no DNA yield | Always dispense exactly the buffer volumes given in the protocol! |
| | • Always follow closely the given instructions with regard to order and mode of mixing (shaking, vortexing, etc). |
| | Add the indicated volume of ethanol (96–100%) to Wash Buffer SW2 Concentrate and mix thoroughly (see section 3 for more information). |
| | Store kit components at room temperature (18-25 °C). Storage at lower temperatures may cause salt precipitation. Check Lysis Buffer SL1 and SL2 for white precipitate. If precipitation occurred, incubate the bottle for 10 min at 30-40 °C and shake every 2 minutes until all precipitate is dissolved (see section 3 for more information). |
| | • Keep bottles tightly closed in order to prevent evaporation or contamination. |
| | Sample material not stored properly |
| | • Whenever possible, use fresh material. |

| Problem | Possible cause and suggestions |
|--|--|
| | Too harsh mechanical sample disruption |
| DNA is degraded | • Reduce intensity or incubation time of mechanical sample lysis. |
| DNA is degraded | DNA is degraded by DNases |
| | • Add at least 10–15 µL Enhancer SX to the lysate. |
| | DNA yield was overestimated |
| | • If DNA eluates are not completely free of contaminants (e.g., RNA, protein, humic substances) UV-VIS quantification based on A_{260} is not reliable due to the contribution of the contaminants to the absorption at 260 nm |
| | Carry-over of ethanol or salt |
| Suboptimal | Make sure to dry the silica membrane and the NucleoSpin[®] Soil Column completely before elution to avoid carry-over of ethanolic Wash Buffer SW2. |
| performance of DNA in downstream | Check if Buffer SW2 has been equilibrated to room temperature (18-25 °C) before use. Washing at lower temperatures decreases the efficiency of salt removal. |
| experiments | Contamination with PCR inhibitors |
| | • The DNA purity can be increased by lowering the amount of starting material (see section 2.4 for more information). |
| | • Enhancer SX can facilitate the release of humic substances. Reduce Enhancer SX to 10 μL or omit the buffer entirely (see section 2.5 for more information). |
| | • Make sure to carefully follow the washing instructions. |
| | • Dilute DNA 1:10 to reduce concentration of inhibitors. |
| | |

6.2 Ordering information

| Product | REF | Pack of |
|---|--------------------|-----------------|
| NucleoSpin [®] Soil | 740780.10/.50/.250 | 10/50/250 preps |
| NucleoSpin [®] DNA Stool | 740472.10/.50 | 10/50 preps |
| NucleoSpin [®] Microbial DNA | 740235.10/.50/.250 | 10/50/250 preps |
| NucleoSpin [®] DNA Lipid Tissue | 740471.10/.50 | 10/50 preps |
| NucleoSpin [®] DNA Insect | 740470.10/.50 | 10/50 preps |
| Buffer SB | 740785.50 | 60 mL |
| Buffer SL1 | 740781.30 | 30 mL |
| Buffer SL2 | 740782.30 | 30 mL |
| Buffer SL3 | 740783.50 | 50 mL |
| Enhancer SX | 740784.50 | 50 mL |
| MN Bead Tube Holder | 740469 | 1 piece |
| MN Bead Tubes Type A (0.6 – 0.8 mm ceramic beads, recommended for soil and sediments) | 740786.50 | 50 pieces |
| MN Bead Tubes Type B (40–400 µm glass beads, recommended for bacteria) | 740812.50 | 50 pieces |
| MN Bead Tubes Type C (1 – 3 mm corundum, recommended for yeast) | 740813.50 | 50 pieces |
| MN Bead Tubes Type D (3 mm steel beads, recommended for insects) | 740814.50 | 50 pieces |
| MN Bead Tubes Type E (40 – 400 µm glass beads and 3 mm steel beads, recommended for hard-to-lyse bacteria within insect samples) | 740815.50 | 50 pieces |
| MN Bead Tubes Type F (1 – 3 mm corundum and 3 mm steel beads, recommended for challenging tissues, e.g., spleen, or lung tissue) | 740816.50 | 50 pieces |
| MN Bead Tubes Type G (5 mm steel beads, recommended for plant material) | 740817.50 | 50 pieces |

| Product | REF | Pack of |
|---|----------------|---------|
| Collection Tubes (2 mL) | 740600 | 1000 |
| MN Bead Tubes 5 mL Type A e.g., 5 mL Bead Tubes containing 0.6–0.8 mm ceramic beads; suitable in conjunction with MN Bead Tube Holder 5 mL or mixer mill* | 740799.50 | 50 |
| MN Bead Tube Holder 5 mL e.g., 5 mL Tube Holder for Vortex-Genie [®] instrument and a 3-inch platform in order to house up to 5 bead tubes | 740459 | 1 |
| NucleoSpin [®] Inhibitor Removal Columns | 740789.50/.250 | 50/250 |

Visit *www.mn-net.com* for more detailed product information.

6.3 Product use restriction/warranty

All MACHEREY-NAGEL products are designed for their intended use only. They are not intended to be used for any other purpose. The description of the intended use of the products can be found in the original MACHEREY-NAGEL product leaflets. Before using our products, please observe the instructions for use and the safety instructions from the respective Material Safety Data Sheet of the product.

This MACHEREY-NAGEL product is carrying documentation stating specifications and other technical information. MACHEREY-NAGEL warrants to meet the stated specifications. The provided warranty is limited to the data specifications and descriptions as given in the original MACHEREY-NAGEL literature. No other statements or representations, written or oral, by MACHEREY-NAGEL's employees, agents or representatives, except written statements signed by a duly authorized officer of MACHEREY-NAGEL are authorized. They should not be relied upon by the costumer and are not a part of a contract of sale or of this warranty.

Liability for all possible damages that occur in any connection with our products is limited to the utmost minimum as stated in the general business terms and conditions of MACHEREY-NAGEL in their latest edition which can be taken from the company's website. MACHEREY-NAGEL does not assume any further warranty.

Products and their application are subject to change. Therefore, please contact our Technical Service Team for the latest information on MACHEREY-NAGEL products. You may also contact your local distributor for general scientific information. Descriptions in MACHEREY-NAGEL literature are provided for informational purposes only.


Last updated: 08/2022, Rev. 04

Please contact: MACHEREY-NAGEL GmbH & Co. KG Tel.: +49 24 21 969-333 support@mn-net.com

Trademarks:

FastPrep is a registered trademark of MP Biomedicals, LLC NucleoSpin[®] is a registered trademark of MACHEREY-NAGEL GmbH & Co KG Vortex-Genie is a registered trademark of Scientific Industries, Inc.

All used names and denotations can be brands, trademarks, or registered labels of their respective owner – also if they are not special denotation. To mention products and brands is only a kind of information (i.e., it does not offend against trademarks and brands and can not be seen as a kind of recommendation or assessment). Regarding these products or services we can not grant any guarantees regarding selection, efficiency, or operation.

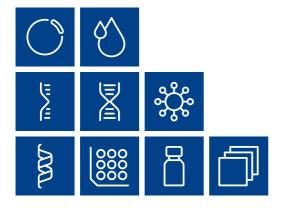

www.mn-net.com

MACHEREY-NAGEL



MACHEREY-NAGEL GmbH & Co. KG DE Tel.: +49 24 21 969-0 info@mn-net.com Valencienner Str. 11 52355 Düren · Germany

CH Tel.: +41 62 388 55 00 sales-ch@mn-net.com FR Tel.: +33 388 68 22 68 sales-fr@mn-net.com US Tel.: +1 888 321 62 24 sales-us@mn-net.com



A038626/0231